- | | |
|-----------|--------------------------------|
| DE | Bedienungsanleitung |
| GB | Instruction manual |
| FR | Mode d'emploi |
| ES | Manual de instrucciones |

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CAUTION!

To work with this microscope, sharp and pointed aids are being used. Please take care that this microscope and its accessories are stored at a place out of reach of children. Let children only work with this microscope under an adult's supervision!

Keep packing material (plastic bags etc.) away from children!

1. Scope of delivery

Stereo microscope (A), interchangeable objectives (B): 1.0x (Art. No. 5802520) or 0.5x/1.0x/2.0x (Art. No. 5802530), 1 pair of WF10x eyepieces (C), 1 pair of rubber eyecups (D), power cable (E), spare fuse (F), dust cover (G), Instruction manual (H)

Also required (not included):

small Philipps screwdriver, small flathead screwdriver

2. Parts overview

- | | |
|--------------------------------------|-------------------------------|
| ① Eyepieces | ⑨ Connecting screw |
| ② Diopter adjustment ring | ⑩ Lock nut |
| ③ Objective with clamping screw (3a) | ⑪ Focusing wheel |
| ④ Front microscope arm | ⑫ Illumination |
| ⑤ Microscope pole | ⑬ Power socket |
| ⑥ Rear microscope arm | ⑭ Fuse holder |
| ⑦ Locking ring | ⑮ On/Off switch |
| ⑧ Microscope base | ⑯ Rubber eyecup |
| | ⑰ replacement fuse (F2AL250V) |

3. Finding the right location for observation

Dry and dust free, room temperature between -5 and +40°C.

Position your device so that it can be disconnected from the power supply at any time. The wall socket should always be located near the device and be easily accessible, since the plug on the power cord serves as a disconnecting device for the power supply.

4. Assembly

1. Place the base (8) on a stable surface.
2. Attach the rear microscope arm (6) to the pole of the microscope base.
3. Use the locking ring (7) to secure the microscope arm.
4. Connect the front part of the microscope arm (4) to the rear arm. Insert the connecting screw (9) through the hole in the rear microscope arm. Take care that the flattened edge shows backwards. Now attach the front microscope arm to the screw. Tighten the lock nut (10) with the lever until the front microscope arm does not tilt down.

5. Connect the power cable to the power connector (13) of the microscope and a free power outlet.

5. Observation

Turn on the microscope using the power switch (15).

Now, look through the eyepieces (1) of the binocular viewing and change the spacing between the eyepieces so that you get a round, clear field of view. The definition is adjusted by slowly turning the focusing wheel (11).

If you have to observe a very large or very small object, you can adjust the microscope. Loosen the holding screw on the side and move appropriately the microscope head on its guidance. If you have found the correct position (distance of 22 cm from the object), tighten the holding screw again.

6. Diopter adjustment

Turn the left diopter adjustment ring (2) on the left eyepiece to the central position (silver marking). Shut your left eye and look with the right eye through the right eyepiece. Now focus with the focusing wheel (11), focus correctly for your right eye. Now, look with your left eye through the left eyepiece. If the image isn't sharp, adjust the focus for your left eye by turning the diopter adjustment ring. Now, the microscope is correctly adjusted.

7. Replacing the objective

Note: The objective of this stereo microscope can be replaced. Depending on the model, additional interchangeable objectives are included or can be purchased as optional accessories.

Hold the mounted objective (3) firmly with one hand and use a small flathead screwdriver to loosen the slotted screw (3a). Remove the objective from the holder. Place the desired interchangeable objective on the holder and screw it on with the slotted screw as well.

8. Replacing the eyepieces

Note: The eyepieces of this stereo microscope can be replaced. Depending on the model, additional eyepieces are included or can be purchased as optional accessories.

Use a small Phillips screwdriver to unscrew the cross screws on the two eyepiece holders. Remove the eyepieces from the holders and insert others. Screw them on with the cross screws as well.

9. Exchange of the fuse

Switch off the microscope and pull out the plug of the power outlet. Grip with a small screwdriver under the lug of the fuse holder (14) and pull it carefully forward. Take out the fuse and change it into a new one (17). Install the fuse holder and plug again.

10. Care and maintenance

10.1. Cleaning the lenses

Sweep the lens with a lens tissue or soft fabric immersed with a small amount of alcohol (e.g. isopropanol).

10.2. Cleaning the painted parts

The dust on the painted parts can be removed by cleaning cloth. For the grease spots, the gauze immersed slightly with aviation gasoline is recommended. Do not use organic solvents such as alcohol, ether or other thinner etc. for cleaning the painted parts or plastic components. Your microscope is a high-quality optical instrument. Avoid disassembling the microscope. Improper disassembling affects the high utility of the microscope. Moreover, the warranty expires. Cover the microscope with the dust protective cover and place it there, where it is dry and dust-free. We suggest the storage of all objectives and eyepieces in a closed container with drying agent.

11. Technical data

	Art. No. 5802520	Art. No. 5802530
Magnification:	10x	5x/10x/20x
Working distance:	230 mm	250 mm (5x)/230 mm (10x)/80 mm (20x)
Power supply:	230V / 50HZ via power cable	
Surge protection:	with fine fuse, type F2AL250V	
Illumination:	LED, long-life, not exchangeable	
Eyepiece holder:	30.5 mm	

12. Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electric equipment together with household waste material! In observance of European Directive 2012/19/EC on waste electrical and electronic equipment (WEEE) and its implementation in accordance with national law, electric equipment that have reached the end of its life must be collected separately and returned to an environmentally compatible recycling facility.

13. Declaration of Conformity

13.1. CE Declaration of Conformity

CE Bresser GmbH has formed a „CE Conformity Declaration“ in line with the applicable directive. This declaration is available for review under https://www.bresser.de/download/5802520/CE/5802520_CE.pdf
https://www.bresser.de/download/5802530/CE/5802530_CE.pdf

13.2. UKCA Declaration of Conformity

**UK
CA** Bresser UK Ltd. has formed a „UKCA Declaration of Conformity“ in line with the applicable directive. This declaration is available for review under https://www.bresser.de/download/5802520/UKCA/5802520_UKCA.pdf
https://www.bresser.de/download/5802530/UKCA/5802530_UKCA.pdf

14. Warranty & Service

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at www.bresser.de/warranty_terms.

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